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English

Note: This section is an English instruction content using the illustration in the first column on the left as an example.

Product Accessories List

- 1.1 Electronic Eyepiece
1.2 Smartphone Stretch Clip
1.3 USB Type-C Charging Cable
1.4 EVA Ring (for connecting and fixing small-sized eyepieces)
1.5 Cleaning Cloth

Introduction to the Electronic Eyepiece

- 2.01 Power and WiFi Switch
2.02 WiFi Indicator Light
2.03 TF Card Slot
2.04 Charging/Working Status Indicator Light
2.05 USB Type C Interface
2.06 Reset Button
2.07 Smartphone Clip Slot
2.08 Smartphone Clip Release Switch
2.09 Lens
2.10 Clip Left and Right Pressure Switches

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Eyepiece Usage Instructions

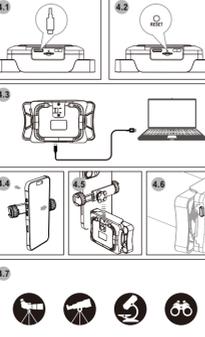
- 3.1 Insert the TF card and press the button for 3 seconds to turn on the device.
3.2 Scan the QR code in the manual with your smartphone to download the APVision software.
3.3 After installation, open the APP and click "Connect Device" to go to the device search WiFi hotspot interface.
3.4 Electronic Eyepiece Indicator Light Description
3.5 When the WiFi connection is successful, it will jump to the real-time preview screen of the APP.
3.6 Connecting to a Telescope
3.7 Connecting to a Microscope
3.8 Connecting to a Microscope

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Other Instructions

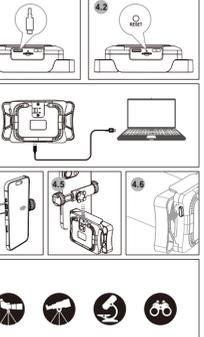
- 4.1 Charging
4.2 Reset
4.3 Data Transfer
4.4 Connecting to a Microscope
4.5 Connecting to a Microscope
4.6 Connecting to a Microscope
4.7 Connecting to a Microscope
4.8 Connecting to a Microscope

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Thank you very much for choosing our smartphone electronic eyepiece. Please read the manual carefully before using this lens and we suggest you keep it well for future reference.

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Please read this manual carefully before choosing our smartphone electronic eyepiece, and we suggest you keep it well for future reference.

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FCC Caution:

This device complies with part 15 of the FCC Rules. Operation is subject to the following two conditions: (1) This device may not cause harmful interference, and (2) this device must accept any interference received, including interference that may cause undesired operation.

Any Changes or modifications not expressly approved by the party responsible for compliance could void the user's authority to operate the equipment.

Note: This equipment has been tested and found to comply with the limits for a Class B digital device, pursuant to part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference in a residential installation. This equipment generates uses and can radiate radio frequency energy and, if not installed and used in accordance with the instructions, may cause harmful interference to radio communications. However, there is no guarantee that interference will not occur in a particular installation. If this equipment does cause harmful interference to radio or television reception, which can be determined by turning the equipment off and on, the user is encouraged to try to correct the interference by one or more of the following measures:

- Reorient or relocate the receiving antenna.
- Increase the separation between the equipment and receiver.
- Connect the equipment into an outlet on a circuit different from that to which the receiver is connected.
- Consult the dealer or an experienced radio/TV technician for help.

The device has been evaluated to meet general RF exposure requirement. The device can be used in portable exposure condition without restriction.