Limited Pre-Series

English

Operating Instructions **Quantos
HPD**

Deutsch

Bedienungsanleitung **Quantos HPD**

Français

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
操作说明书 **Quantos HPD**

METTLER TOLEDO

The logo graphic consists of a series of parallel, slightly curved lines that form a diamond shape, with the text 'METTLER TOLEDO' centered over it.

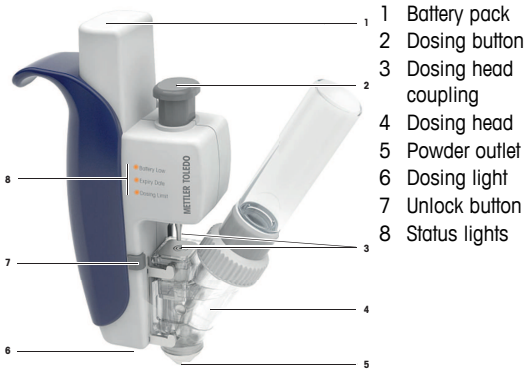1 Overview

2 Safety Information

2.1 Definition of Warnings and Symbols

General hazard

CAUTION

for a hazardous situation with low risk, resulting in damage to the device or the property or in loss of data, or minor or medium injuries if not avoided.

2.2 Product Specific Safety Notes

CAUTION

Damage of Device

- For use only in dry interior rooms.
- Use only the original batteries in the battery pack delivered with your instrument.
- Do not open the instrument: It does not contain any parts which can be maintained, repaired, or replaced by the user. If you ever have problems with your instrument, contact your METTLER TOLEDO dealer.

Always operate and use your instrument only in accordance with the instructions contained in this manual. The instructions for setting up your new instrument must be strictly observed.

If the instrument is not used according to these Operating Instructions, protection of the instrument may be impaired and METTLER TOLEDO assumes no liability.

Intended Use

Your Dosing System is used for weighing and dosing powder or liquid samples. Use the instrument exclusively for this purpose. Any other type of use and operation beyond the limits of technical specifications without written consent from Mettler-Toledo AG, is considered as not intended.

Certain samples require special care!

With certain types of samples, there is a possibility of danger to personnel or damage of property. Please note that the user always has the responsibility and liability for damage caused by use of any types of samples!

3 Operation

3.1 Switching on and off

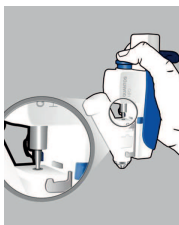
The instrument switches on automatically as soon as you pick it up. The instruments can also be switched on by pressing the dosing button.

The instrument will automatically switch off if you put down the instrument for more than 15 seconds or if you do not press the dosing button for 45 seconds.

3.2 Dosing

- Insert dosing head.
⇒ Dosing light turns on.

- To couple, press once the dosing button.

- Press the dosing button beyond coupling position.
⇒ Powder outlet opens.
⇒ Stirring and tapping will be actuated and powder flows.

If the dosing head which you insert is read and found to be valid for use then the dosing light (white light next to the Powder Open-

ing) will light faintly. As soon as the motor begins to turn then the light intensity increases and illuminates the opening of the vial into which you want to dose.

Releasing coupling

- To release coupling, press the small unlock button once.

Releasing dosing head

- To release dosing head, press the small unlock button during removing the dosing head.

3.3 Status lights

The status lights are three orange LEDs. They can only be seen when lit which indicate the following criteria:

LED-Name	Blinking (Warning)	Permanent (No more doses possible)
Battery Low	approx. 10% life time remaining	Empty
Expiry Date	Will be reached within next 5 days	Expired zero
Dosing Limit	5 doses remaining	Zero doses remaining

Blinking LED when the motor is blocked

If you press the dosing button but the motor does not start and the dosing light does not shine, then the corresponding LED will blink rapidly, because one of the above listed criteria is invalid. When you release the dosing button the LED will be permanently lit.

Expiry Date

For each dosing head you can set the **Expiry Date** of the powder. To set this you need a Quantos system with RFID reading/writing unit. If no expiry date is set then the corresponding LED will remain inactive. If you have set an expiry date and the dosing head date has expired then you will not be able to perform further doses.

Dosing Limit

The **Dosing Limit** is like a dosing counter. Each dosing head has a limited number of doses the so called dosing limit. The dosing limit can be set by the user within pre-set boundaries. The factory setting dosing heads is 250 doses (from a range from 1 to 999 doses). For the sample dosing head the factory setting is 36 doses (from a range of 1 to 99 doses). To change the factory settings you need a Quantos system with RFID reading/writing unit.

4 Maintenance

4.1 Cleaning

CAUTION

Damage of instrument

- Ensure that NO liquid comes in contact with the battery pack or the dosing unit.
- Read the instruction on the battery pack before inserting the batteries.
- Use the delivered batteries.
- Do NOT mix different types or brands of batteries.

Your instrument is made from high quality, resistant materials and can therefore be cleaned with a commercially available, mild cleaning agent and damp cloth.

4.2 Changing batteries

Replace batteries in the battery pack when indicated with the status light **Battery Low**.

Order number 30066628

- 1 Remove battery pack.
- 2 Remove old batteries and insert new batteries described on the battery pack. (one additional pair of batteries is delivered with this package).
- 3 Install battery pack. Check that there is NO gap between battery pack and instrument.

4.3 Disposal

In conformance with the European Directive 2002/96/EC on Waste Electrical and Electronic Equipment (WEEE) this device may not be disposed of in domestic waste. This also applies to countries outside the EU, per their specific requirements.

Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. If you have any questions, please contact the responsible authority or the distributor from which you purchased this device. Should this device be passed on to other parties (for private or professional use), the content of this regulation must also be related.

Thank you for your contribution to environmental protection.

5 Technical Data


General Data

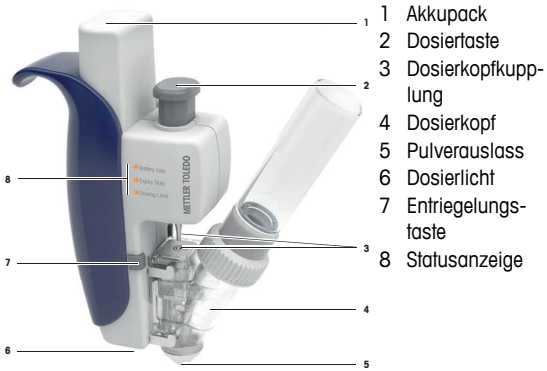Power Supply:	Battery pack with 6 V Lithium batteries, containing 2 pieces CR123A, 3 V, 1400 mAh
Shortcut Protection:	PTC High Current Discharge Protection: 5 A
Net weight:	184 g (without dosing head)
Degree of pollution:	Level 2
Standards for safety and EMC:	See Declaration of Conformity
Range of application:	For use only in enclosed, dry interior rooms

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© Mettler-Toledo AG 03/2015
Mettler-Toledo AG, Laboratory Weighing